## Oligo (dT)<sub>15</sub> Primer

### **Product Description:**

A primer designed to initiate synthesis of a cDNA from total RNA in a reverse transcription reaction. For use in generating labeled cDNA for screening microarrays.

**Form:** Supplied in lyophilized form

**Storage:** 4°C or -20°C

**MW:** 4501 mol

**T<sub>m</sub>:** 26°C

**Purity:** Purified by reverse phase chromatography, >95%

**Pack size:** 30 µg equals ~ 1 OD unit (A260). 30 µg -6.7 nmol

**Quality Control:** Purity is verified by analytical PAGE

### **Notes for Use**

Stored at 4°C or at -20°C when lyophilised; once the product is resuspended in nuclease-free water store at **-20°C**. For maximum recovery prior removing the cap of the tube spin it briefly to bring down the content. Reconstitute with nuclease-free water (included vial) and vortex briefly. Avoid repeat freeze-thaw cycles. Aliquots are stable for at least 24 months at -20°C. Further dilutions can be made in assay buffer.

### **Order Information:**

| Ref.   | Product                  |       |
|--------|--------------------------|-------|
| 20.401 | Random Primers           | 30 µg |
| 20.402 | Oligo (dT) <sub>15</sub> | 30 µg |

### **Notice to buyers/users:**

Some of the applications which may be performed with this product are covered by applicable patents in certain countries. The purchase of this product does not include or provide a license to perform patented applications. Users may be required to obtain a license depending on the country and/or application.

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