

---

# **Termi-DNA-Tor** - Complete removal of bacteria, mycobacteria, fungi, yeast, virus and nucleic acids (DNA or RNA) from working surfaces

**STORE AT 25°C**

---

## ***DESCRIPTION***

Regular cleaning and disinfection of laboratory working surfaces from contaminants is a must in molecular biology, and laboratories working with microorganisms. The choice of active ingredient is very important because determine its exposure time and effectiveness. The current employed methods are time consuming and toxic, moreover its long term use lead to a deterioration of cleaned surfaces.

## ***PROPERTIES***

**Termi-DNA-Tor** makes a difference in everyday cleaning routine, as it is an efficient tool for complete **removal of bacteria, mycobacteria, fungi, yeast, virus and nucleic acids (DNA or RNA)** present in the laboratory environment. Its regular use keeps working surfaces free of contaminants, thus minimising risk of false results, while minimally affecting the cleaned surfaces<sup>1</sup>.

Use of Termi-DNA-Tor is direct and easy, as its aerosol presentation allows achievement of the best level of cleanliness while using the minimum amount of product.

## ***INSTRUCTIONS FOR USE***

Ready-to-use solution. Professional use

Wash previously the working surface with a cleaning agent, spray the solution directly over the surface to treat, until it's completely wet and let it act for the time required for the desired antimicrobial activity

EFFECTIVENESS	BACTERICIDE	MYCOBACTERICIDE	YEASTICIDAL	FUNGICIDE	ACTIVE AGAINST VIRUS
<b>Exposure Time</b>	30 Seconds	5 Minutes	30 Seconds	5 Minutes	30 Seconds

---

<sup>1</sup> For surfaces sensitive to ethanol try Termi-DNA-Tor first in a small area of the surface

## **FORMULATION**

Hydroalcoholic solution (55% etanol), quaternary ammonium propionate, perfume

## **SAFETY INSTRUCTIONS**

- Very flammable liquid and steam. It causes severe eye irritation.
- Avoid contact with skin and eyes.
- Wear suitable gloves.
- In case of contact with eyes, rinse immediately with plenty of water for several minutes and seek medical advice.
- Keep away from food, drinks and animal feeding stuffs.
- Avoid use on materials sensitive to alcohol.
- Do not use near sparks and open flames
- Do not spray on hot surfaces.
- In case of accident or if you feel unwell, seek medical advice immediately (show this document if possible).
- Keep out of children's reach..
- Do not light flames or smoke during application.
- Store it in a well ventilated place
- Avoid breathing spray mist
- Harmful to aquatic organisms causing long-term adverse effects. Avoid release to the environment
- Don't reuse packaging
- This product and its packaging are classified as hazardous waste
- Store at (5°C – 25°C)
- More information available on the MSDS sheet of the product



## **ORDER INFORMATION**

<b>Cat. No.</b>	<b>Product</b>
22.001	Termi-DNA-Tor (450 ml)
22.002	Termi-DNA-Tor (2 x 450 ml)

**EXCLUSIVELY FOR PROFESSIONAL USE**